

Ansell

HOW TO CONNECT A GLOVE TO AN ISOLATOR SLEEVE - INSTRUCTION SHEET

Please use the instructions below as a guide to attaching a glove (e.g BFAP) to an isolator sleeve (e.g. GSL33NITPP26) with a BBCO-100 connector.



1 Prepare the isolator sleeve and connector ready for assembly, with the sleeve inside out



2 Take the sleeve and slide the connector over the bottom of it




3 Now, pull the bottom of the sleeve over the connector and tuck under the connector so that the sleeve is secured in place



4 Take the glove and feed it through the cuff connector, fingers first



5 Pull the cuff of the glove over the connector

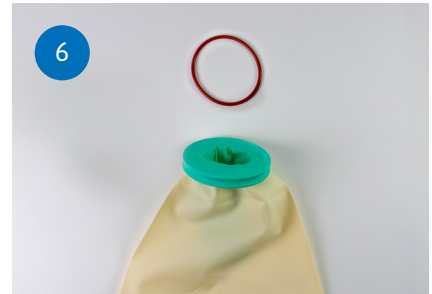

6 Continue as in photo #5, ensuring to completely cover the sleeve cuff connector with the cuff of the glove



7 Take the silicone rubber O-ring and fix into groove of connector over the glove



8 Turn the assembled sleeve and glove the correct way round



9 Finally, place hand into fully assembled sleeve/glove system